




The MICROSCOOP[®] MINT System Grant Support Document

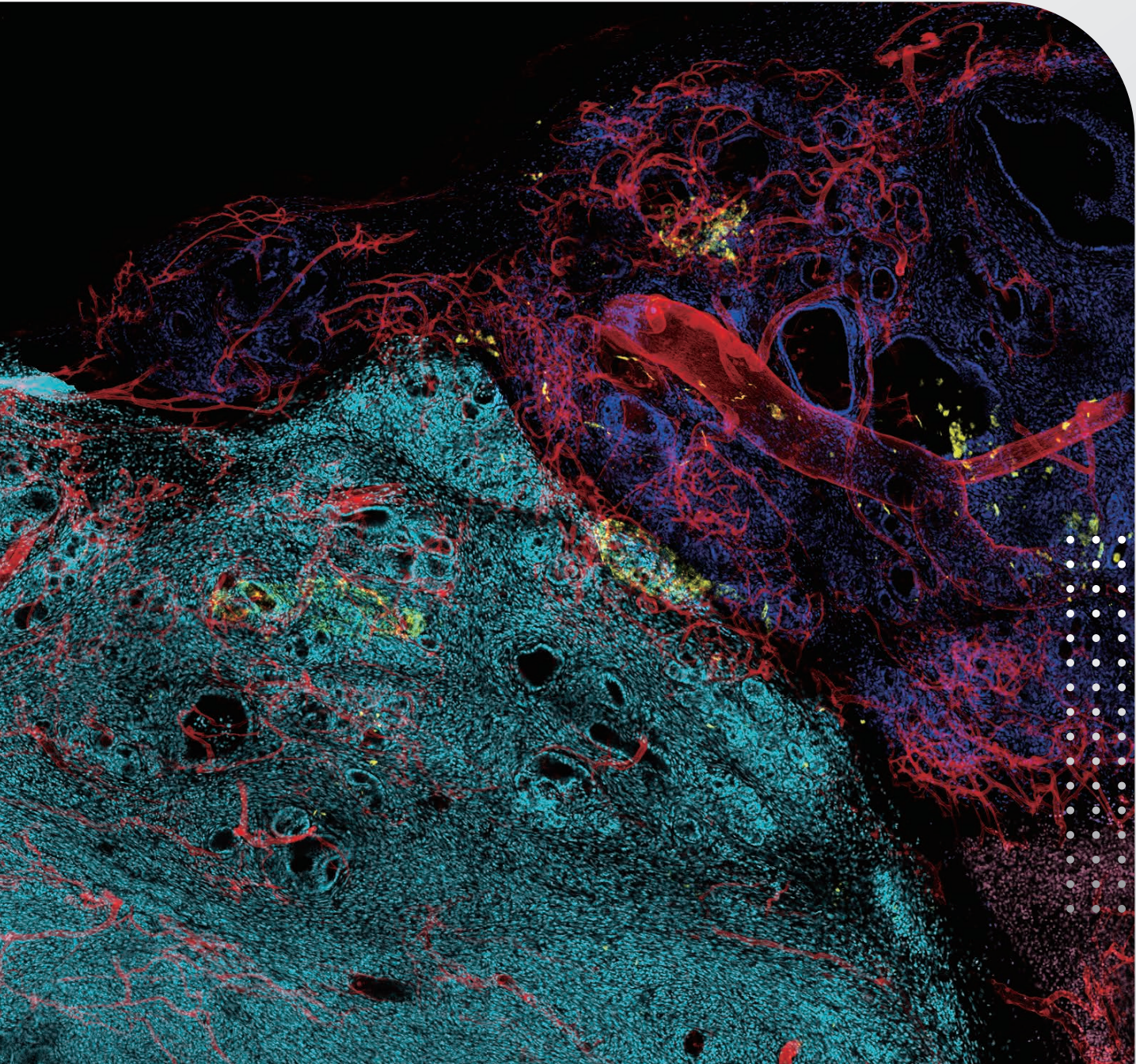


Syncell Inc. approves the use of images, graphs/charts, methods, and descriptions found in this grant package with proper citing for grant proposal purposes. Please cite: [Image, data, methods, description] courtesy of Syncell Inc.

Research use only. Not for use in diagnostic procedure.

Why Spatial Discovery?

Spatial discovery represents a significant shift in proteomic research, focusing on the de novo identification of proteins in specific cell compartments and microenvironments. Unlike spatial targeted profiling that rely on predefined sets of known proteins that require antibodies or aptamers, spatial discovery starts with using Microscoop® to precisely extract the proteins localized to selected regions of cells, cellular compartments, or structures. These extracted proteins are then subjected to mass spectrometry analysis for identification and characterization.

The key advantage of spatial discovery lies in its ability to unveil previously unrecognized proteins that may play crucial roles in cellular processes within specific spatial contexts. By casting a wide net for protein identification, spatial discovery offers a comprehensive view of the protein landscape within selected cellular regions at subcellular precision, opening new avenues for understanding cellular function and signaling pathways.

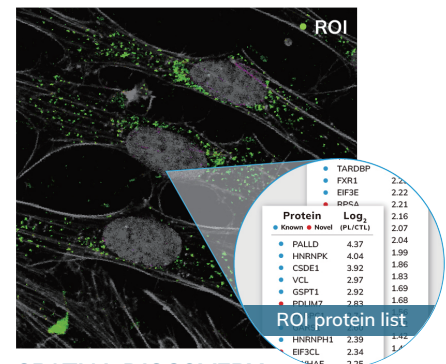

SPATIAL DISCOVERY
Identifying proteins in regions of interest (ROIs)

Platform Overview

Microscoop® Mint is the solution for spatial protein purification.


- Isolates proteins from specific subcellular locations with high precision.
- Enables unbiased discovery of new protein constituents at your region of interest (ROI).
- ROI's can be generated at the level of hundreds of nanometers.
- No protein candidate list is needed in advance.
- Automatically photo-biotinylates proteins in millions of similar spatial targets.
- Discovers novel biomarkers or therapeutic targets of disease-associated locations.
- Subsequent LC-MS/MS analysis to reveal novel spatial proteomes.



How Microscop[®] Works


STEP1 Real-Time Image Analysis

Photolabeling kit (i.e. Synlight-Pure[™] Kit or Synlight-Rich[™] Kit) is first added to a cell or tissue (fixed, FFPE or frozen) sample for a photochemical reaction. An image mask is created from a user-defined stain (IF, brightfield, and more) for regions of interest (ROI's) that define the local area of interest for where proteins are extracted. After the sample is loaded onto the stage, Microscop[®] takes an image (or images of multiple colors) of the sample at one field of view (FOV) at a time. The image or images are analyzed in real time by Microscop's software Autoscoop[™], which executes traditional image processing or AI deep learning to segment the user's region of interest. Pre- or post-processing can be included to enhance segmentation accuracy.

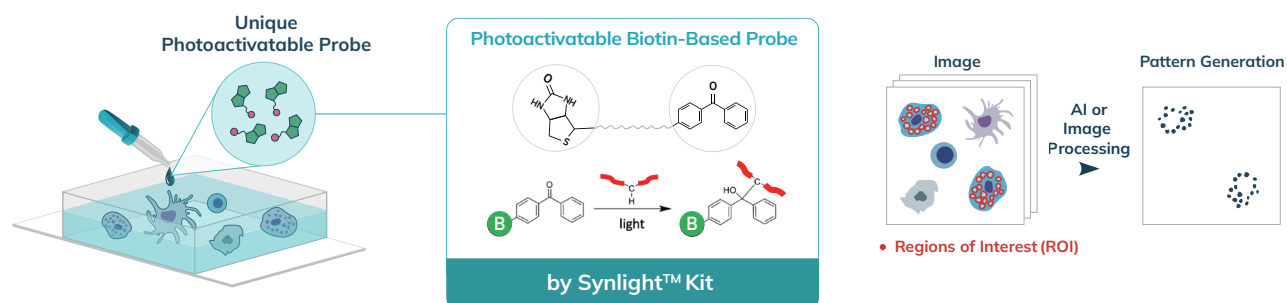


STEP2 Patterned Photo-Biotinylation

A femtosecond light source is controlled to illuminate the segmented region of interest one point at a time. This patterned illumination triggers targeted protein photo-biotinylation in high spatial precision through the reactions of light-sensitive probes of Synlight-Pure[™] Kit or Synlight-Rich[™] Kit. This patterned photolabeling is repeated for thousands of FOVs automatically to assure that enough proteins are biotinylated for later proteomics analysis using mass spectrometry.

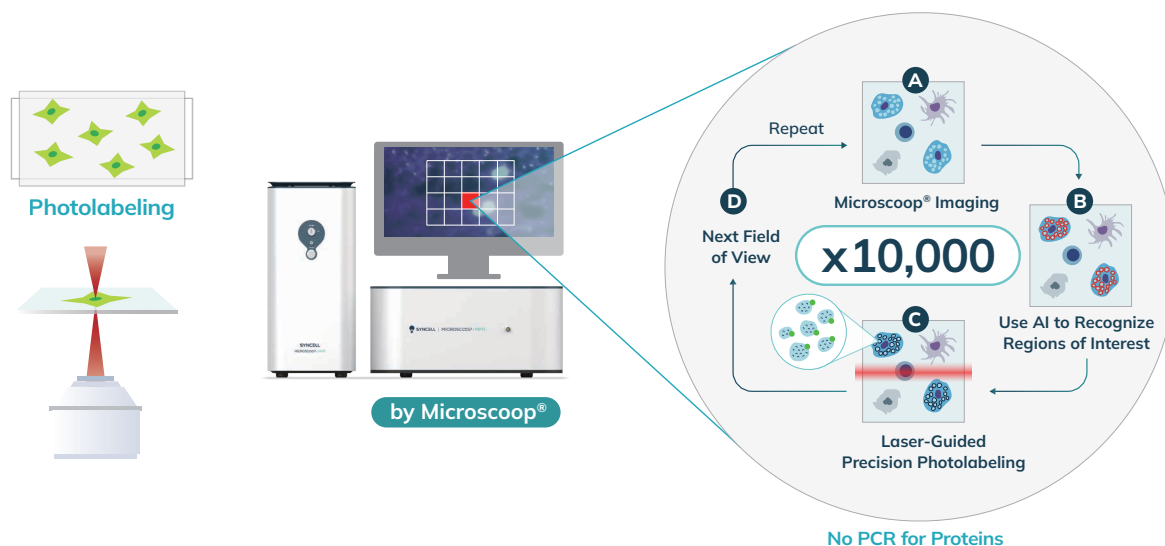


STEP3 Protein Extraction

Photolabeled cells or tissues are scraped from the slide or chamber. Materials from multiple slides or chambers can be pooled together to increase the total protein contents. The scraped sample is then treated with reagents of protein extraction kit (i.e. Synpull™) Kit to lyse the sample, enrich the proteins by immunoprecipitation, and digest them into peptides for proteomics analysis.

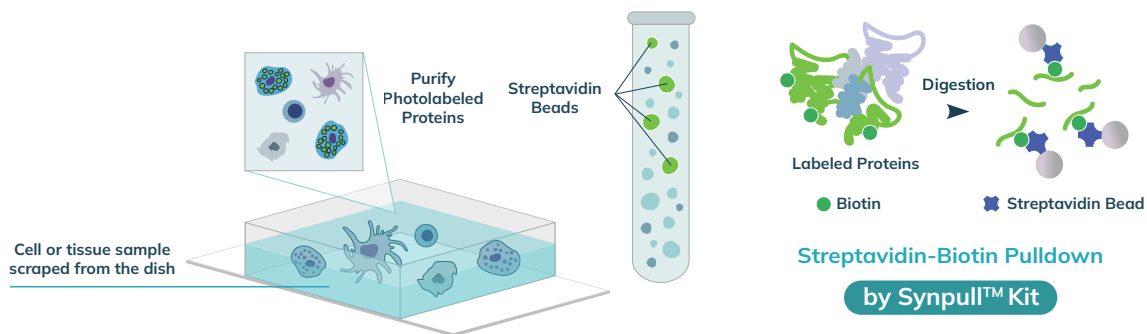

STEP4 Proteomic Identification

The collected peptides are sent to a mass spectrometer to perform LC-MS/MS analysis. Proteomes of both the photo-labeled and unlabeled (CTL) samples are obtained. By comparing the control and photolabeled proteomes, a location-specific proteome at the region of interest is obtained in high sensitivity, high specificity, and high spatial precision. Validation can be done by colocalization of immunostaining or additional functional assays.

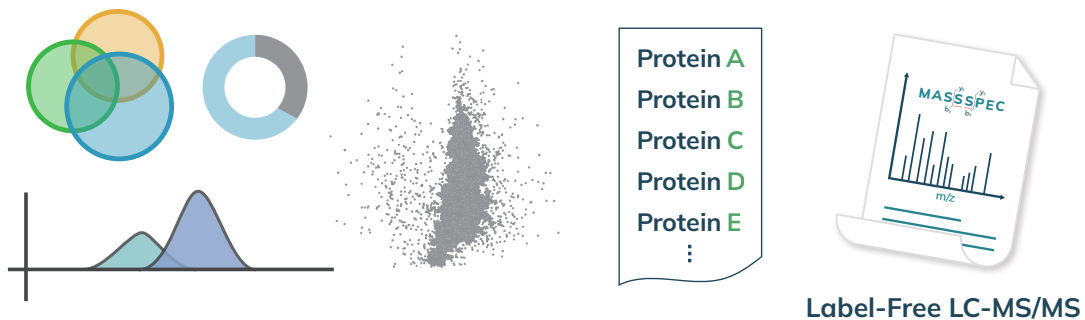

The Microscope[®] Mint Specifications

DESCRIPTION		SPECIFICATIONS
Function		Optoproteomics: Ultra-content high-speed microscopy-guided subcellular photoaffinity labeling for hypothesis-free high-precision proteomic discovery
Workflow		Cyclic procedure of the following step: 1. Microscopy imaging: image acquisition 2. Pattern segmentation: selection of user-defined regions of interest 3. Patterned scanning illumination: point-by-point photochemical reactions 4. Stage movement: change of the field of view
Components		<ol style="list-style-type: none"> 1 Microscope[®] system (optical engine and electrical controller) 2 Inverted epifluorescence microscope Filter sets for microscope 3 Epifluorescence illumination light source 4 Two-photon laser for Microscope[®] photolabeling 5 Camera 6 Software package
1	Dimensions (L x W x H)	Electrical controller : 44 cm x 22 cm x 47 cm Optical engine: 68 cm x 46 cm x 22 cm
	Power Source	100 - 240 VAC, 50/60 Hz
2	Objectives	10x (up to NA 0.45) 20x (up to NA 0.80) 40x (up to NA 0.95)
	Stage	Motorized XY positioning stage (X: ±57 mm, Y: ±36.5 mm stroke) with a vessel holder, suitable for microscope slides, chamber slides, or micro-dishes
3	Imagery Wavelength	Dyes: e.g. DAPI, FITC, Cy3, Cy5 Fluorescent proteins: e.g. EBFP2, EGFP, DsRed/mCherry
5	Camera	sCMOS camera (resolution: 2048 x 2048, pixel size: 6.5 µm x 6.5 µm)
	Binning Options	Low resolution mode: 800 x 800 pixels High resolution mode: 1600 x 1600 pixels
6	Operating System	Microsoft Windows 10
	Pattern Segmentation Options	Toolbox for traditional image processing Trained model using AI deep learning
Labeling Resolution		300 nm+**
Sample Format		Cells - fixed on a chambered coverslip Tissues - slide mounted FFPE (5 - 10 µm in thickness) or frozen tissue section (10 - 20 µm in thickness)
Sample Size Requirement		Cell numbers: 4 x 10 ⁵ to 1 x 10 ⁶ cells for a single LC-MS/MS analysis* Tissue slides: 4 - 8 tissue section for a single LC-MS/MS analysis*

*Application dependent on and varying with the area and the number of ROIs. **Objective dependent

All product specifications and data are subject to change without notice to improve reliability, function, design, or otherwise. For Research Use Only. Not for use in diagnostic procedures.

Supporting Data

Reproducibility of Microscoop®

SynCell Microscoop® demonstrates a robust identification and analysis of stress granules (SGs). Accurate photolabeling and efficient image processing techniques were developed, resulting in high reproducibility of SG image masking, photo-labeling, and proteomics experiments. Novel SG proteins were identified and validated, contributing to a better understanding of SG composition.

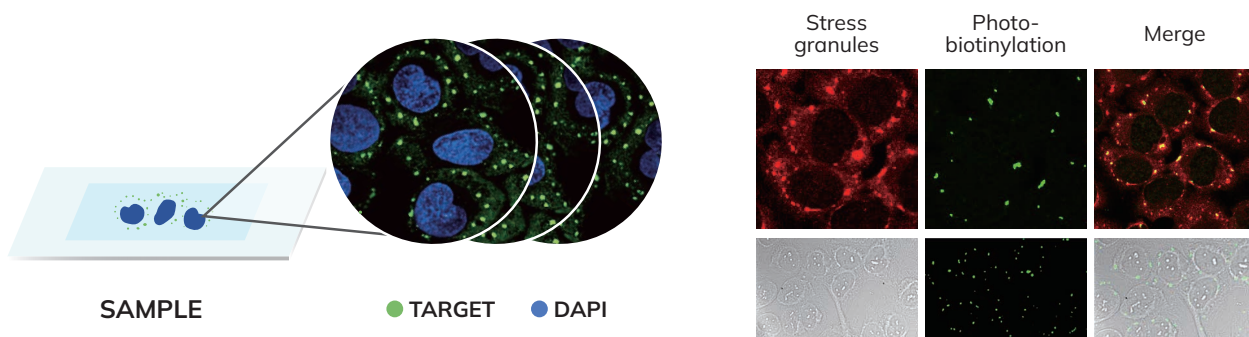


FIGURE 1 | Precise and accurate photolabeling for each stress granule in U-2OS cells.

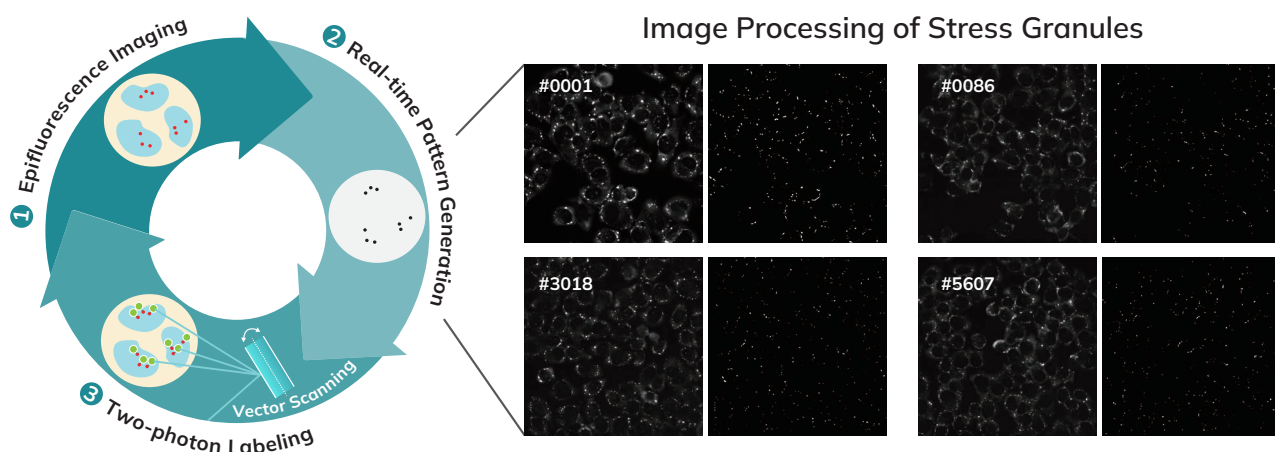


FIGURE 2 | Robust image processing isolates stress granules from background noise in each field of view (FOV), generating consistent patterns in all FOVs.


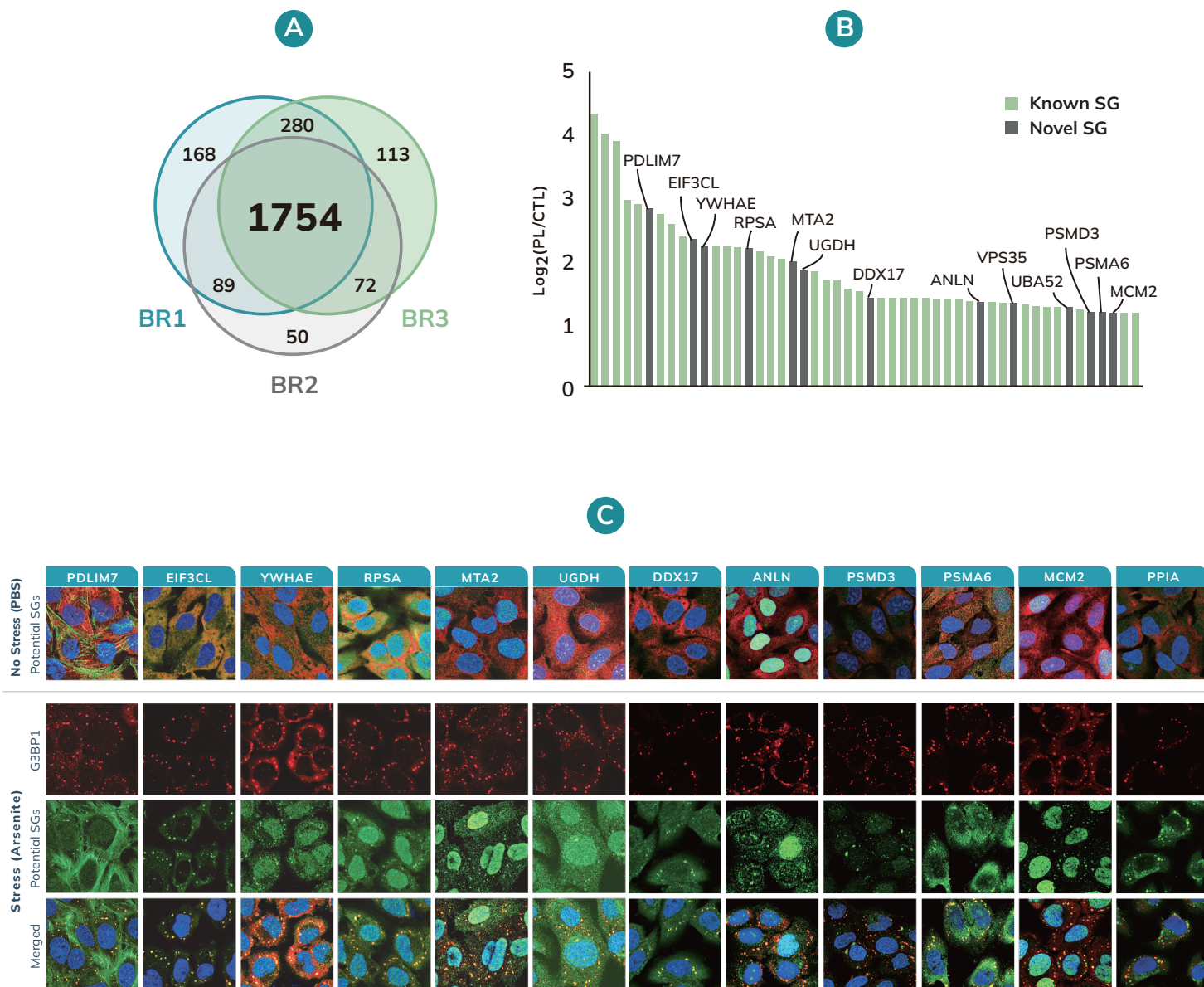


FIGURE 3 | A, High reproducibility was demonstrated between three biological replicates of spatial photolabeling SG proteomics. B, Identification of potentially novel SG protein among top-ranked spatially purified proteins. C, Validation of novel SG proteins by immunostaining in U-2OS cells, which form SGs under arsenite stress. Proteins identified as potential SG proteins (green) are highly co-localized with SG marker G3BP1 (red).

Applications

Microscope® is a microscopy-guided protein-picking system designed for precise photobiotinylation at up to 240nm resolution. If you can clearly define your region of interest under a microscope using immunofluorescent imaging, Microscope® can achieve targeted protein labeling. It supports a variety of dyes, including DAPI, FITC, Cy3, Cy5, and fluorescent proteins like EBFP2, EGFP, and DsRed/mCherry. Traditional image processing tools, such as adaptive thresholding and filtering, can be employed to define specific regions of interest. Applicable cellular patterns include subcellular structures, cell-cell contact sites, organelle contact sites, entire cells, and aggregates. Microscope® has been successfully used with formalin-fixed, paraffin-embedded (FFPE), fresh frozen tissue samples, and fixed cultured and primary cell samples.

Broad Discover Applications

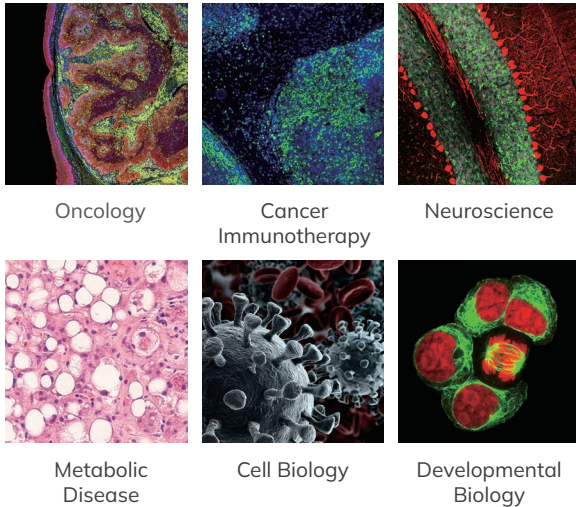

Oncology

Cancer Immunotherapy


Neuroscience

Metabolic Disease

Cell Biology

Developmental Biology

Applicable Cellular Patterns

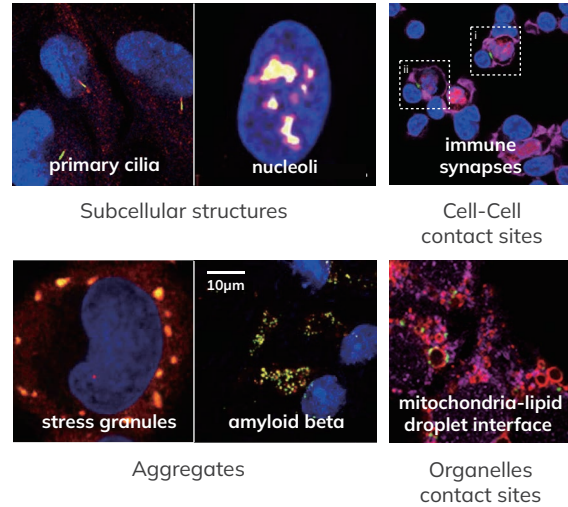

primary cilia

nucleoli

immune synapses

Subcellular structures

Cell-Cell contact sites

stress granules


amyloid beta

mitochondria-lipid droplet interface

Aggregates

Organelles contact sites

Applicable Sample Types

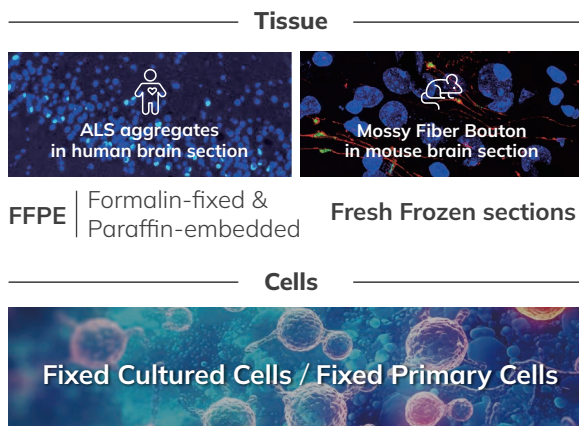

Tissue

ALS aggregates in human brain section

Mossy Fiber Bouton in mouse brain section

FFPE

Formalin-fixed & Paraffin-embedded

Fresh Frozen sections

Cells

Fixed Cultured Cells / Fixed Primary Cells

Applicable Areas



Disease Mechanisms
Novel protein discovery

Microscope® Consumables

SynCell offers optimized reagent kits for Microscope®. Synlight Kits are the proprietary kits that contain photochemical probes and other needed reagents for photolabeling. Synlight-Rich™ Kit is suitable for features larger than 300 nm with high labeling efficiency. Synpull™ Kit contains a large group of reagents optimized for low-volume pulldown and mass spectrometry-ready preparation suitable for both cell and tissue samples.


The following content is commercially available for precise photolabeling and pulldown:

Photolabeling Kit

Synlight-Rich™ Kit


Up to 6 reactions (1-3 rounds** of LC-MS/MS)

*Positive control included **1 round=1 test group+1 control group

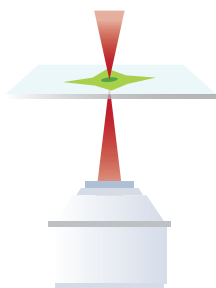

Targeted photolabeling by
MICROSCOOP®

Synlight-Rich™ enables high efficiency photolabeling, suitable for sub-micron or micron size structures, such as the cell bodies (~10 μm) and axons (~2 μm) of Purkinje cells as well as nucleoli (~1 μm).

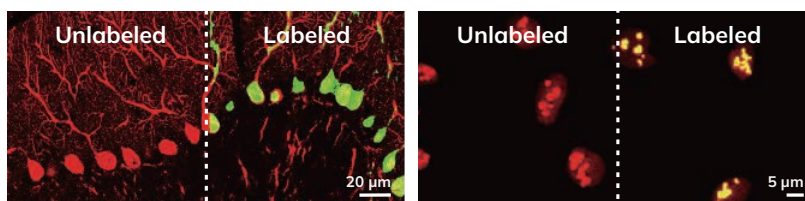

Purkinje cells
in the mouse cerebellum


Nucleoli

Protein Extraction Kit


Synpull™ Kit

Up to 6 reactions (3 rounds** of LC-MS/MS)

**1 round=1 test group+1 control group

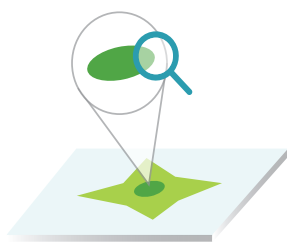

Extraction and processing
of photolabeled samples
for MS-based proteomics

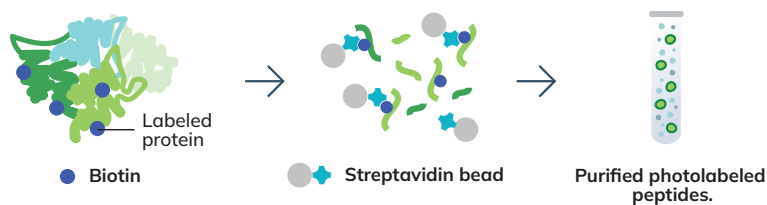

Benefits to consider for your research

High Sensitivity

Microscoop® mechatronics controls the illumination of photoreactive reagents to accurately label all proteins in user specified locations.

High-Content

Users can perform multi-omics spatial labeling to identify and analyze tens and thousands of individual cells.

High Specificity

High Specificity-Unique combination of computer vision imaging and accurate photolabeling provides unsurpassed high specificity for protein biomarker discovery from a particular cellular or subcellular structure.

AI-Driven

Custom machine learning software solutions allow users to define different cellular and subcellular structures.

Photolabeling

Microscoop® utilizes two-photon excitation microscopy to provide accurate XYZ position labeling with subcellular resolution.

Speed

Imaging takes 0.1 seconds, followed by AI or image processing, which requires 0.1 to 0.5 seconds since the calculations are performed in real time.

Additional Resources

Training

The Microscoop® Mint support team offers expert guidance throughout the entire user experience, from installation to ongoing training. Users are invited to participate in virtual training sessions that cover a range of topics.

Additionally, Microscoop® users benefit from exclusive access to user-group meetings and workshops, where they can exchange best practices and success stories within the community, fostering a collaborative environment for maximizing the technology's potential.

Literature

For more details on Microscoop® platform, refer to our publications and supporting literature.

[Publication](#)

[Product literature and research posters](#)

Additional Support

General inquiries: info@syncell.com

Learn more about spatial protein purification solutions: [contact a sales representative](#)

Phone: +1 617-631-2746

