



SYNCELL GRASP Cell Sample Preparation and Shipping Guideline

Guideline for cell sample preparation

- Grow cell cultures in 2-well chamber slides.(Do you need to use adherent cells or you can spin down suspension cells?)
- Thickness of chamber slide should less than 170µm
- Need to describe the confluency you need.
- Cells should be fixed by 4% paraformaldehyde (PFA) for 10 minutes.
- **Samples should be immersed in PBS to avoid drying.**

Shipping Considerations

- Pack the slides in a slide box ad seal the outside with parafilm.
- Label the outside of the slide box with following information.
 - Principal Investigator name
 - Email
 - Date
- Place a copy of sample manifest and the filled Customer Information form (see below) in the package.
- Ship the package at ambient temperature.

Customer Information

Name:

Email:

Institution/ Company:

Principal Investigator:

Country:

Address:

Date: